

Vol. EN09



NMR
SHIGEMI®

NMR/EPR Tubes
&
Accessories

SHIGEMI Co., LTD.

2-4-15 Nakanosanno Hachioji
TOKYO JAPAN

Pone: +81-42-624-2207 FAX: +81-42-622-0937 info@shigemi.co.jp
<https://www.shigemi.co.jp/en-introduction>

Contents

1. Symmetrical MICRO NMR Tubes	2~6
2. Thin wall Symmetrical MICRO NMR Tubes	7~10
3. SLOT TUBE	11~12
4. Ultra Precision NMR Tubes(Thin wall)	13
5. Economy NMR Tubes	14
6. 5mm MICRO Bottom NMR Tubes	15
7. 5mm ϕ Coaxial NMR tube set	16~17
8. 5mm ϕ Coaxial INNER TUBE for Heat sealing	18
9. 5mm ϕ Coaxial GAS tight NMR tube set	19~20
10. Quartz NMR Tubes	21
11. Ultra Thin wall NMR Tubes	22~23
12. Precision Large Volume NMR Tubes	24
13. ESR (EPR) Quartz Tubes	25
14. TRUE GAS TIGHT TUBE FOR NMR · ESR	26~27
15. Accessories	28
16. 5mm NMR SAMPLE TUBE for SampleJet	29






Symmetrical MICRO NMR Tubes

These unique NMR microtubes are made out of a special type of hard glass with excellent chemical durability. They are magnetic susceptibility matched to each of the following solvents: CDCL₃, CD₃OD, D₂O, DMSO


Therefore, the best resolution of a sample can be obtained in each of these solutions with minimal amounts of sample. Yet, the use of a small amount of solution will not affect the spectral resolution or the sideband.

How to use Symmetrical MICRO NMR Tubes

- ★ Do not spinning
- ★ We do not garanty auto shimming
Plaese make shim file from 0
- ★ Plaese take out bables
 - ① You turn in an inner tubeslowly
 - ② Brief stop when a inner tube tach at sample surface.
 - ③ You push quick a inner tube
 - ④ Bables was disapred return an inner tube right possition

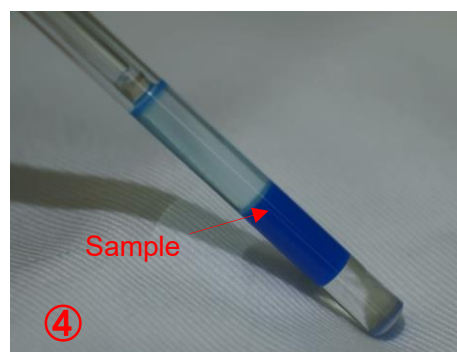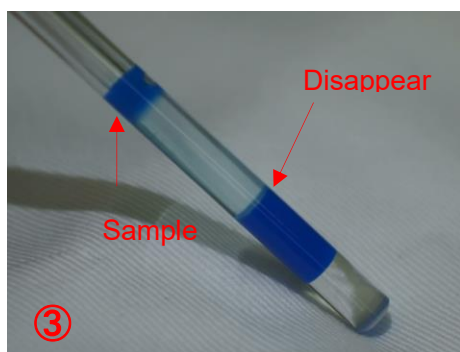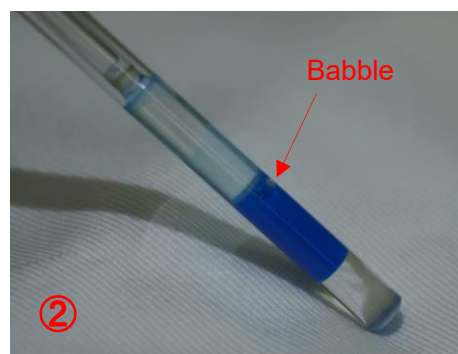


Symmetrical MICRO NMR Tubes for D₂O

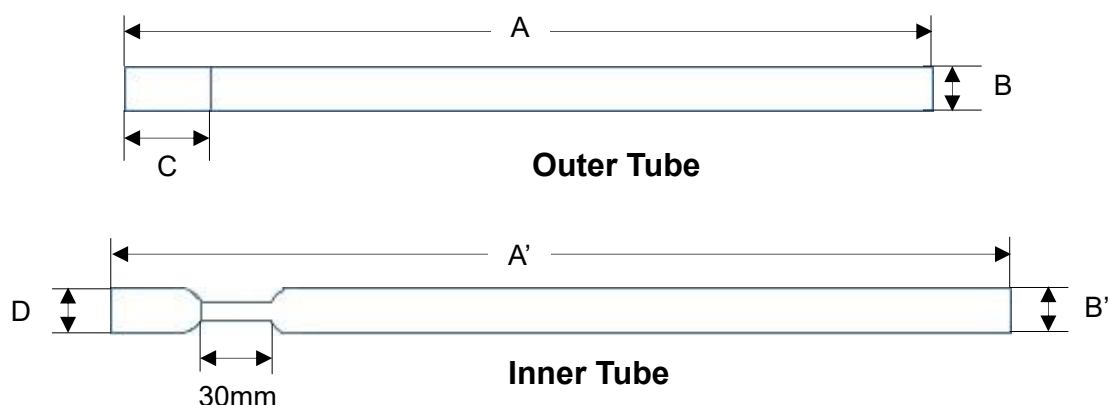

Parts N.O.		Outer Tube(mm)			Inner Tube(mm)			Price 1set = 1 outer tube and 1 inner tube JPY
		Length A	O.D B	Muched Glass C	Length A'	O.D. B'	Muched glass D	
2.5mmφ	BMS-0025	120	2.5 +0 -0.01	6	115	1.8	1.86 +0 -0.01	8,100
3mmφ	BMS-003	180	3.00 -0.01 -0.02	10	190	2.4	2.47 +0 -0.01	9,200
4mmφ	BMS-004B	180	4.000 +0.005 -0.010	8	190	3.1	3.16 +0 -0.01	10,400
	BMS-004J	180	4.000 +0.005 -0.010	12	190	3.1	3.16 +0 -0.01	10,400
	BMS-005V	180	4.000 +0.005 -0.010	15	190	3.1	3.16 +0 -0.01	10,400
5mmφ	BMS-005B	180	4.965 +0.005 -0.010	8	190	4.1	4.17 +0 -0.01	10,400
	BMS-005J	180	4.965 +0.005 -0.010	12	190	4.1	4.17 +0 -0.01	10,400
	BMS-005V	180	4.965 +0.005 -0.010	15	190	4.1	4.17 +0 -0.01	10,400
8mmφ	BMS-008BJ	190	7.98 ±0.01	10	200	6.9	6.97 +0 -0.01	20,700
	BMS-008V	190	7.98 ±0.01	15	200	6.9	6.97 +0 -0.01	23,000
10mmφ	BMS-010BJ	190	9.98 +0 -0.01	10	200	8.9	8.97 +0 -0.01	25,300
	BMS-010V	190	9.98 ±0.01	15	200	8.9	8.97 +0 -0.01	28,800


Symmetrical MICRO NMR Tubes for CDCl₃

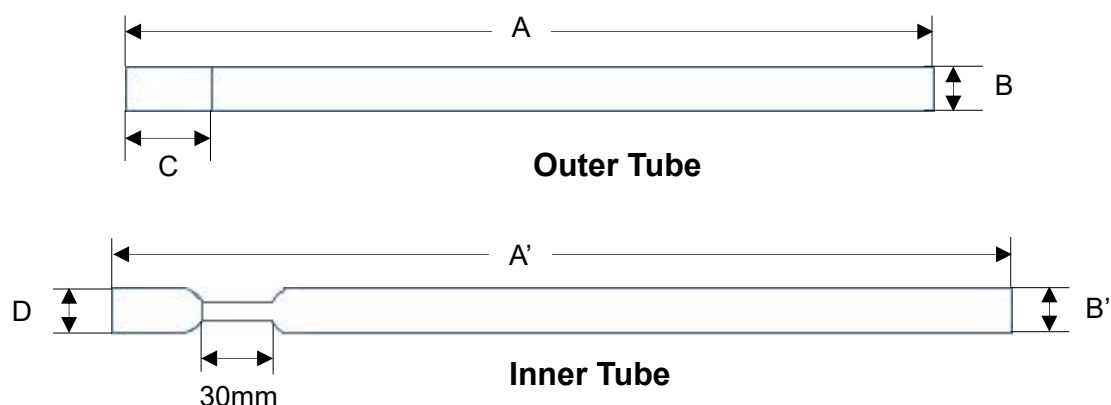

Parts N.O.		Outer Tube(mm)			Inner Tube(mm)			Price 1set = 1 outer tube and 1 inner tube JPY
		Length A	O.D. B	Muched glass C	Length A'	O.D. B'	Muched glass D	
2.5mmφ	CMS-0025	120	2.5 +0 -0.01	6	115	1.8	1.86 +0 -0.01	8,100
3mmφ	CMS-003	180	3.00 -0.01 -0.02	10	190	2.4	2.47 +0 -0.01	9,200
5mmφ	CMS-005B	180	4.965 +0.005 -0.010	8	190	4.1	4.17 +0 -0.01	10,400
	CMS-005J	180	4.965 +0.005 -0.010	12	190	4.1	4.17 +0 -0.01	10,400
	CMS-005V	180	4.965 +0.005 -0.010	15	190	4.1	4.17 +0 -0.01	10,400
8mmφ	CMS-008BJ	190	7.98 ±0.01	10	200	6.9	6.97 +0 -0.01	20,700
	CMS-008V	190	7.98 ±0.01	15	200	6.9	6.97 +0 -0.01	23,000
10mmφ	CMS-010BJ	190	9.98 +0 -0.01	10	200	8.9	8.97 +0 -0.01	25,300
	CMS-010V	190	9.98 ±0.01	15	200	8.9	8.97 +0 -0.01	28,800




Symmetrical MICRO NMR Tubes for DMSO (CD₃)₂SO, C₆D₆, C₅D₅N

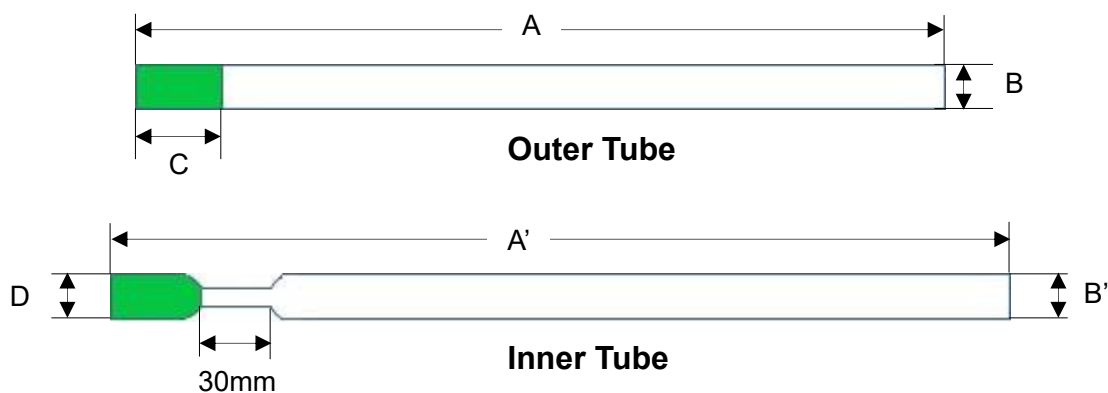

Parts N.O.		Outer Tube(mm)			Inner Tube(mm)			Price 1set = 1 outer tube and 1 inner tube JPY
		Length A	O.D. B	Muched glass C	Length A'	O.D. B'	Muched glass D	
2.5mmφ	DMS-0025	120	2.5 +0 -0.01	6	115	1.8	1.86 +0 -0.01	8,100
3mmφ	DMS-003	180	3.00 -0.01 -0.02	10	190	2.4	2.47 +0 -0.01	9,200
5mmφ	DMS-005B	180	4.965 +0.005 -0.010	8	190	4.1	4.17 +0 -0.01	11,500
	DMS-005J	180	4.965 +0.005 -0.010	12	190	4.1	4.17 +0 -0.01	11,500
	DMS-005V	180	4.965 +0.005 -0.010	15	190	4.1	4.17 +0 -0.01	11,500
8mmφ	DMS-008BJ	190	7.98 ±0.01	10	200	6.9	6.97 +0 -0.01	20,700
	DMS-008V	190	7.98 ±0.01	15	200	6.9	6.97 +0 -0.01	23,000
10mmφ	DMS-010BJ	190	9.98 +0 -0.01	10	200	8.9	8.97 +0 -0.01	25,300
	DMS-010V	190	9.98 ±0.01	15	200	8.9	8.97 +0 -0.01	28,800


Symmetrical MICRO NMR Tubes for CD₃OD, CD₃CN

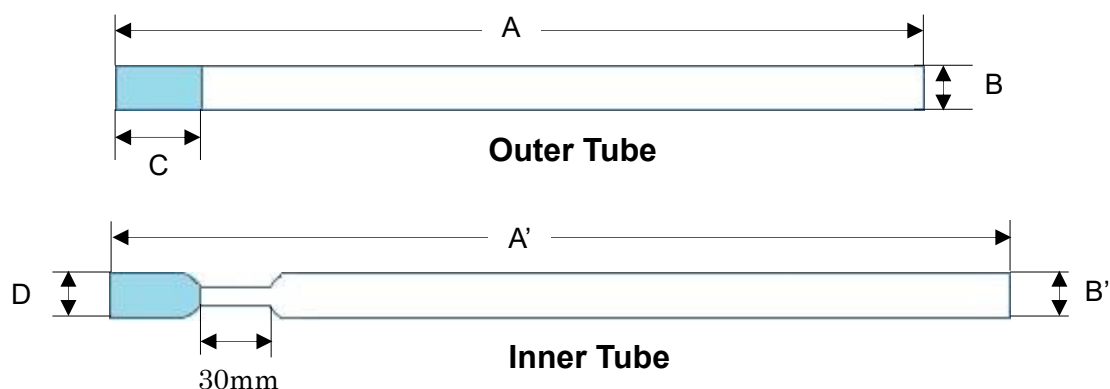

Parts N.O.		Outer Tube(mm)			Inner Tube(mm)			Price 1set = 1 outer tube and 1 inner tube JPY
		Length A	O.D B	Muched glass C	Length A'	O.D. B'	Muched glass D	
2.5mmφ	MMS-0025	120	2.5 +0 -0.01	6	115	1.8	1.86 +0 -0.01	8,100
3mmφ	MMS-003	180	3.00 -0.01 -0.02	10	190	2.4	2.47 +0 -0.01	9,200
5mmφ	MMS-005B	180	4.965 +0.005 -0.010	8	190	4.1	4.17 +0 -0.01	11,500
	MMS-005J	180	4.965 +0.005 -0.010	12	190	4.1	4.17 +0 -0.01	11,500
	MMS-005V	180	4.965 +0.005 -0.010	15	190	4.1	4.17 +0 -0.01	11,500
8mmφ	MMS-008BJ	190	7.98 ±0.01	10	200	6.9	6.97 +0 -0.01	20,700
	MMS-008V	190	7.98 ±0.01	15	200	6.9	6.97 +0 -0.01	23,000
10mmφ	MMS-010BJ	190	9.98 +0 -0.01	10	200	8.9	8.97 +0 -0.01	25,300
	MMS-010V	190	9.98 ±0.01	15	200	8.9	8.97 +0 -0.01	28,800




Thin wall Symmetrical MICRO NMR Tubes for D₂O

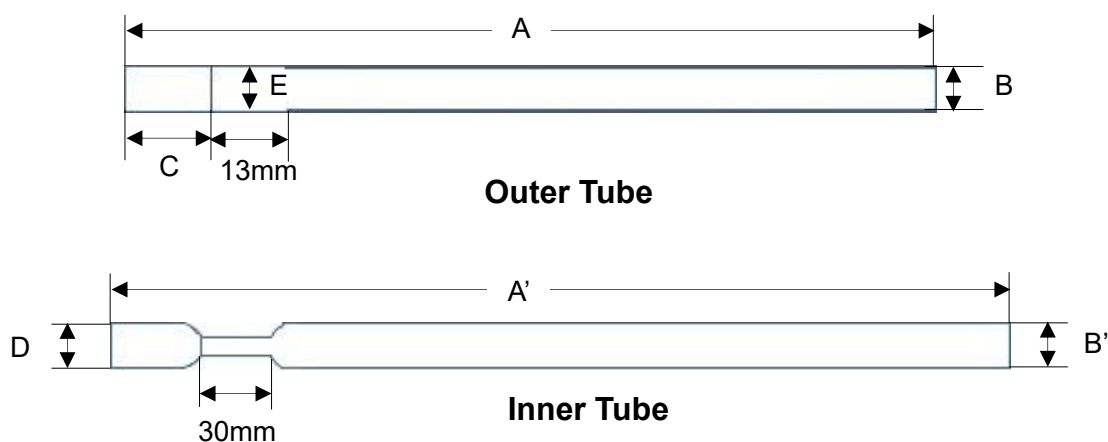


Parts N.O.		Outer Tube(mm)				Inner Tube(mm)			Price(JPY) 1set = 1 outer tube and 1 inner tube
		Length A	O.D. B	Muched Glass C	Thin Wall I.D. E	Length A'	O.D. B'	Mached Glass O.D. D	
5mmφ	BMS-005TB	180	4.965 +0.005 -0.010	8	4.52	190	4.1	4.17 +0 -0.01	11,500
	BMS-005TJ	180	4.965 +0.005 -0.010	12	4.52	190	4.1	4.17 +0 -0.01	11,500
	BMS-005TV	180	4.965 +0.005 -0.010	15	4.52	190	4.1	4.17 +0 -0.01	11,500

Thin wall Symmetrical MICRO NMR Tubes for CDCl_3

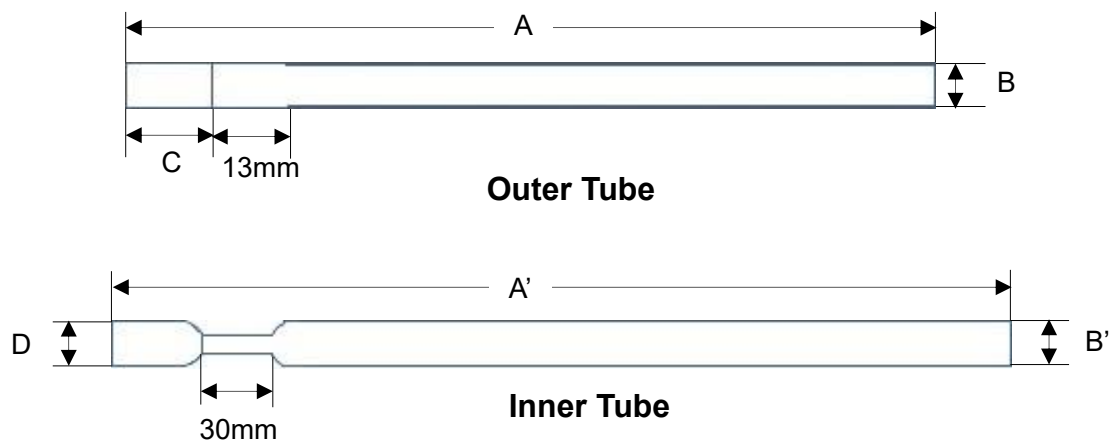

Parts N.O.		Outer Tube(mm)				Inner Tube(mm)			Price(JPY) 1set = 1 outer tube and 1 inner tube
		Length A	O.D. B	Mached Glass C	Thin Wall I.D. E	Length A'	O.D. B'	Mached Glass O.D. D	
5mmφ	CMS-005TB	180	4.965 +0.005 -0.010	8	4.52	190	4.1	4.17 +0 -0.01	11,500
	CMS-005TJ	180	4.965 +0.005 -0.010	12	4.52	190	4.1	4.17 +0 -0.01	11,500
	CMS-005TV	180	4.965 +0.005 -0.010	15	4.52	190	4.1	4.17 +0 -0.01	11,500


Thin wall Symmetrical MICRO NMR Tubes for DMSO (CD₃)₂SO , Benzene C₆D₆ , Pyridine C₅D₅N

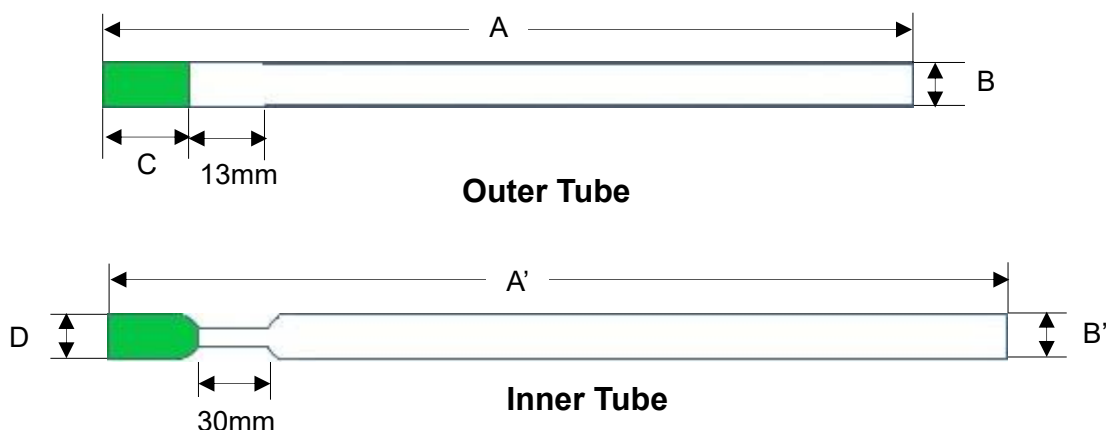

Parts N.O.		Outer Tube(mm)				Inner Tube(mm)			Price(JPY) 1set = 1 outer tube and 1 inner tube
		Length A	O.D. B	Muched Glass C	Thin Wall I.D. E	Length A'	O.D. B'	Mached Glass O.D. D	
5mmφ	DMS-005TB	180	4.965 +0.005 -0.010	8	4.52	190	4.1	4.17 +0 -0.01	12,700
	DMS-005TJ	180	4.965 +0.005 -0.010	12	4.52	190	4.1	4.17 +0 -0.01	12,700
	DMS-005TV	180	4.965 +0.005 -0.010	15	4.52	190	4.1	4.17 +0 -0.01	12,700



Thin wall Symmetrical MICRO NMR Tubes for Methanol CD₃OD, Acetonitrile CD₃CN

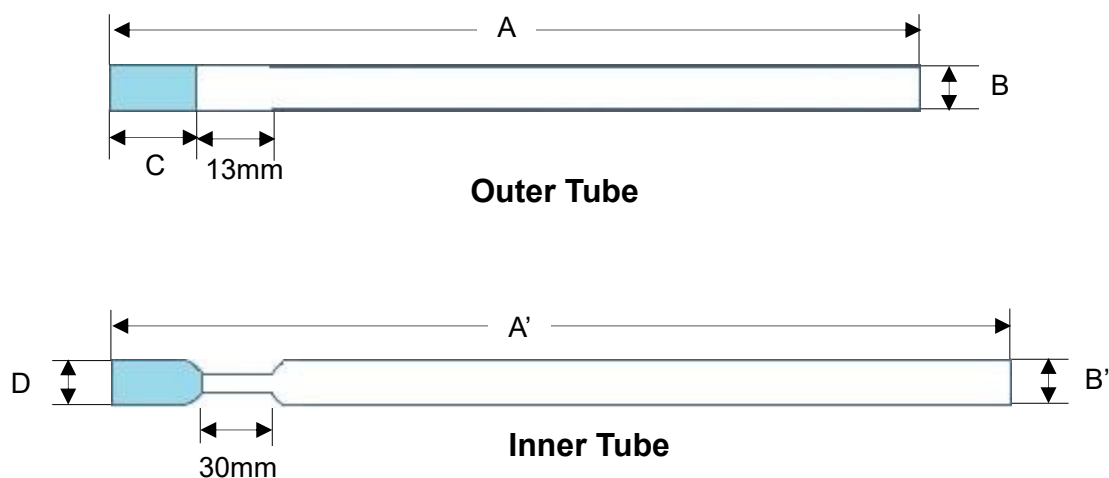

Parts N.O.		Outer Tube(mm)				Inner Tube(mm)			Price(JPY) 1set = 1 outer tube and 1 inner tube
		Length A	O.D. B	Muched Glass C	Thin Wall I.D. E	Length A'	O.D. B'	Mached Glass O.D. D	
5mmφ	MMS-005TB	180	4.965 +0.005 -0.010	8	4.52	190	4.1	4.17 +0 -0.01	12,700
	MMS-005TJ	180	4.965 +0.005 -0.010	12	4.52	190	4.1	4.17 +0 -0.01	12,700
	MMS-005TV	180	4.965 +0.005 -0.010	15	4.52	190	4.1	4.17 +0 -0.01	12,700




SLOT TUBE


For High concentration electrolyte samples

The SHIGEMI Slot Tube provides the highest sensitivity and shortest 90 degree pulse width with any probes when working with lossy samples such as proteins in aqueous solutions containing high concentrations of electrolyte as salt.


Benefit

- ★ Avoid the hot spot of RF power dissipation.
- ★ Able to obtain the best S/N ratio than any other NMR sample tubes.
- ★ Provide the shortest 90 degree pulse width than any NMR sample tubes.
- ★ Has already been confirmed working well with a CryoTM probe (Bruker) 600, 950MHz and a Cold Probe(Varian/Agilent)600MHz.

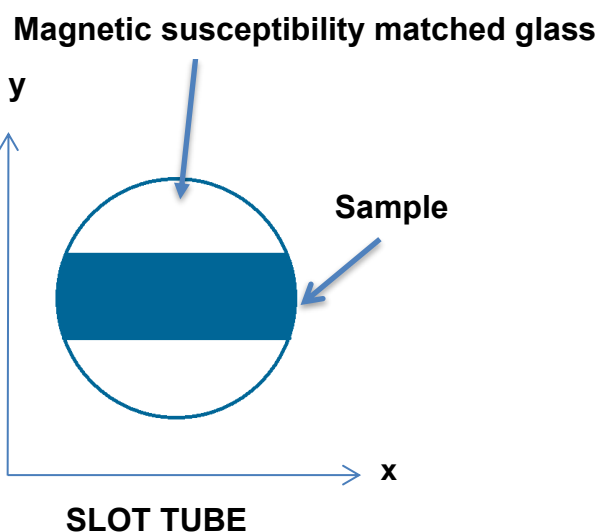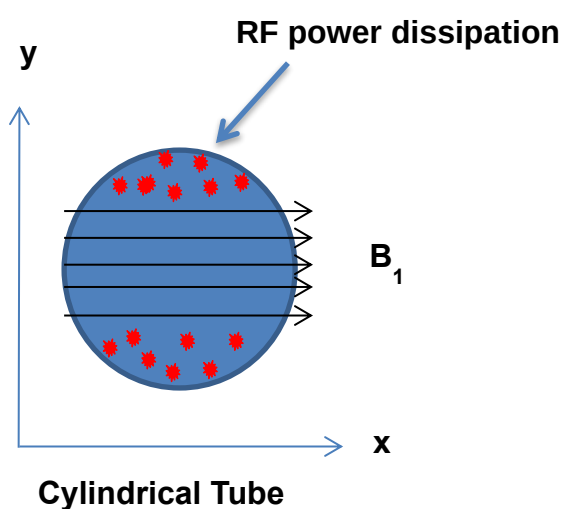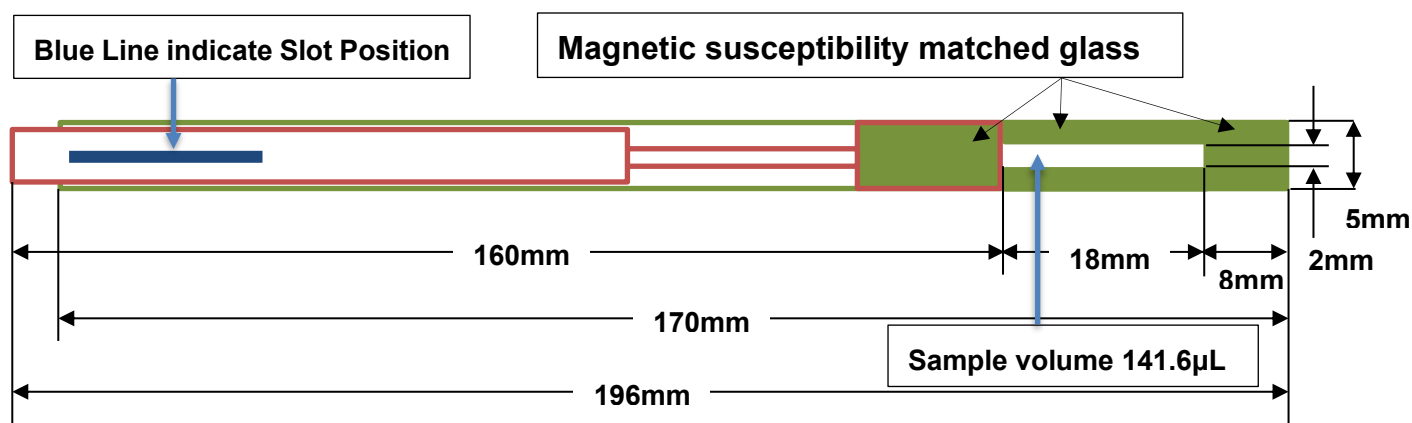

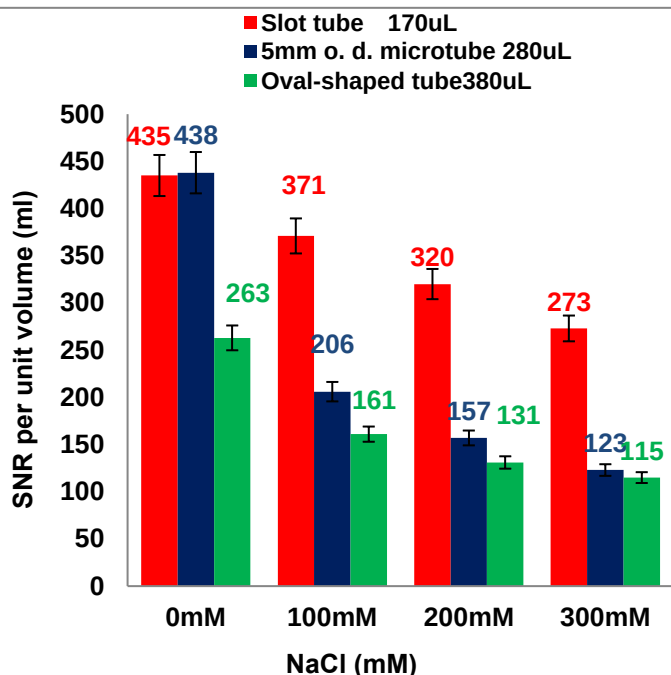

SNR per unit volume(ml) in different tubes
And salt concentrations

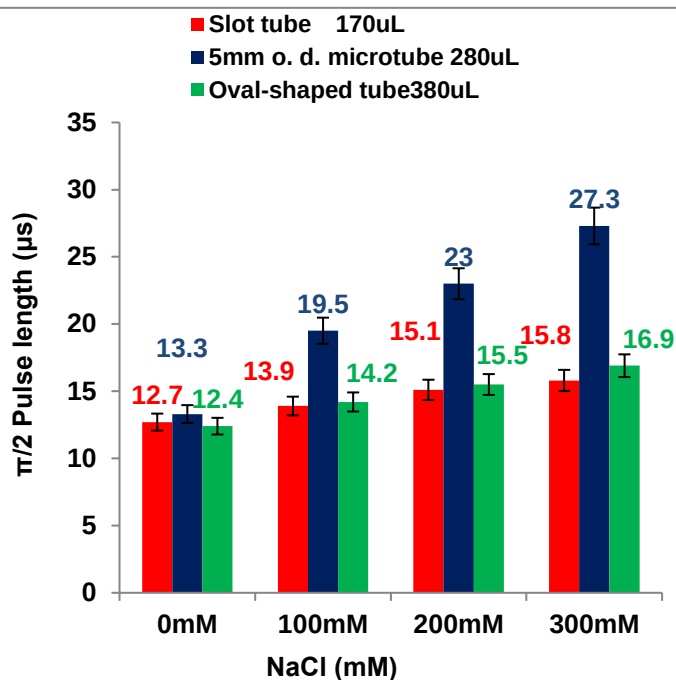

$\pi/2$ Pulse length (μ s) in different tubes
and salt concentrations

Data acquired at 14.1T (600 MHz CryoTMProbe) for the ubiquitin samples of 0.3mM of [¹³C/¹⁵N] containing 10mM sodium phosphate

SLOT TUBE

Magnetic susceptibility matched glass




Slot for sample
144.6μL

SBMS 1 set ¥28,800JPY

Positioning tool

If you do not have “Bruker Shaped Shigemi Tube Shuttle (5mm)”, you need this tool for fixing exact position. (shortest 90 degree puls width)

Do not use the air-lift function.

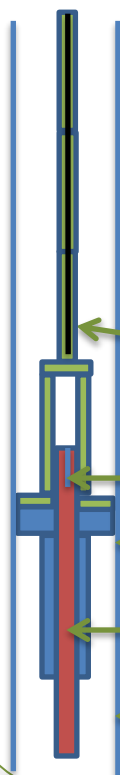

Black Line adjust with Blue Line

Blue Line


Spinner

SLOT TUBE

Upper Barrel

Positioning tool 1 set 172,500JPY

Ultra Precision NMR Sample Tubes

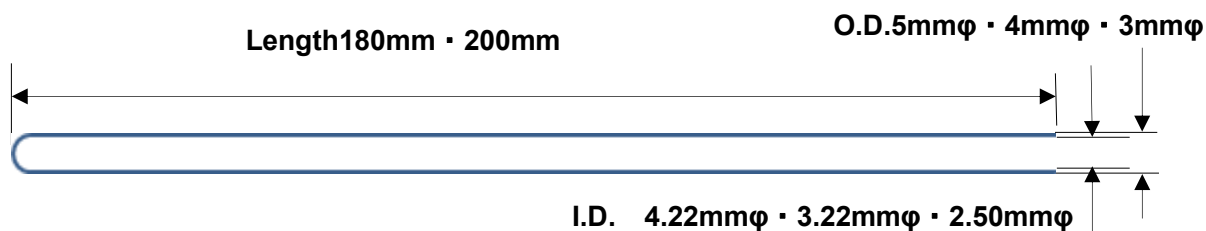

5mmφ Ultra Precision NMR Sample Tubes **Thin wall**

DURAN®

Parts N.O.	Dimension (mm)			Tolerance (μm)		Price (1piece) JPY
	Length	O.D.	I.D.	Straightness	wall thickness control	
PS-001-7	180	4.965 +0.005 -0.010	4.220 +0.01 -0	10	5	2,200
PS-001-8	200	4.965 +0.005 -0.010	4.220 +0.01 -0	10	5	2,300
PS-002-7	180	4.965 +0.005 -0.010	4.220 +0.01 -0	15	10	1,900
PS-002-8	200	4.965 +0.005 -0.010	4.220 +0.01 -0	15	10	2,000
PS-003-7	180	4.965 +0.005 -0.010	4.220 +0.01 -0	15	20	1,700
PS-003-8	200	4.965 +0.005 -0.010	4.220 +0.01 -0	15	20	1,800
PS-004-7	180	4.965 +0.005 -0.010	4.220 +0.01 -0	20	25	1,100
PS-004-8	200	4.965 +0.005 -0.010	4.220 +0.01 -0	20	25	1,200

4mmφ Ultra Precision NMR Sample Tubes **Thin wall**

Parts N.O.	Dimension (mm)			Tolerance (μm)		Price (1piece) JPY
	Length	O.D.	I.D.	Straightness	wall thickness control	
PR-001	180	4 +0 -0.01	3.22 +0.010 -0	15	10	2,300
PR-002	180	4 +0 -0.01	3.22 +0.010 -0	20	20	1,800


3mmφ Ultra Precision NMR Sample Tubes **Thin wall**

Parts N.O.	Dimension (mm)			Tolerance (μm)		Price (1piece) JPY
	Length	O.D.	I.D.	Straightness	wall thickness control	
PN-001	180	3.0 +0.01 -0.02	2.5 +0.01 -0	15	5	2,300
PN-002	180	3.0 +0.01 -0.02	2.5 +0.01 -0	25	20	1,800




Economy NMR Sample Tubes

For routine examination. A quality as same PS-004-7 as Ultra Precision Tubes
We can sale 100 tubes only.

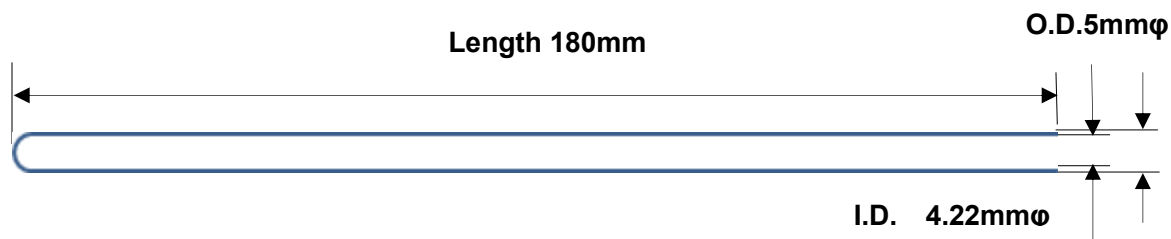

5mm Economy NMR Sample Tubes **Thin Wall** DURAN®

Parts N.O.	Dimension (mm)			Tolerance (μm)		Price JPY	A unit
	Length	O.D.	I.D.	Straightness	Wall thickness control		
PS-005	180	4.965 +0.005 -0.010	4.220 +0.01 -0	25	40	75,000	100 peace

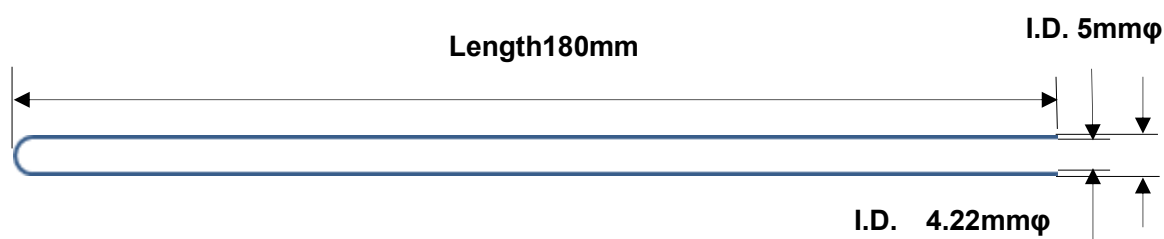

★Matreal is Brosilicate glass ASTM Type1 Class A as DURAN® of Shott AG

★You can use V.T. examination

★Minimum sales unit 1 unit(25pcsX8box)


DURAN®

5mm Disposable NMR Sample Tubes **Thin Wall**

Parts N.O.	Dimension (mm)			Straightness(μm)	price
	Length	O.D.	I.D.		200PCS
EC-57	180	4.965 +0.005 -0.010	4.220 ±0.10	30	44,000
EC-58	200	4.965 +0.005 -0.010	4.220 ±0.10	30	46,000



5mm ϕ Micro Bottom Tubes

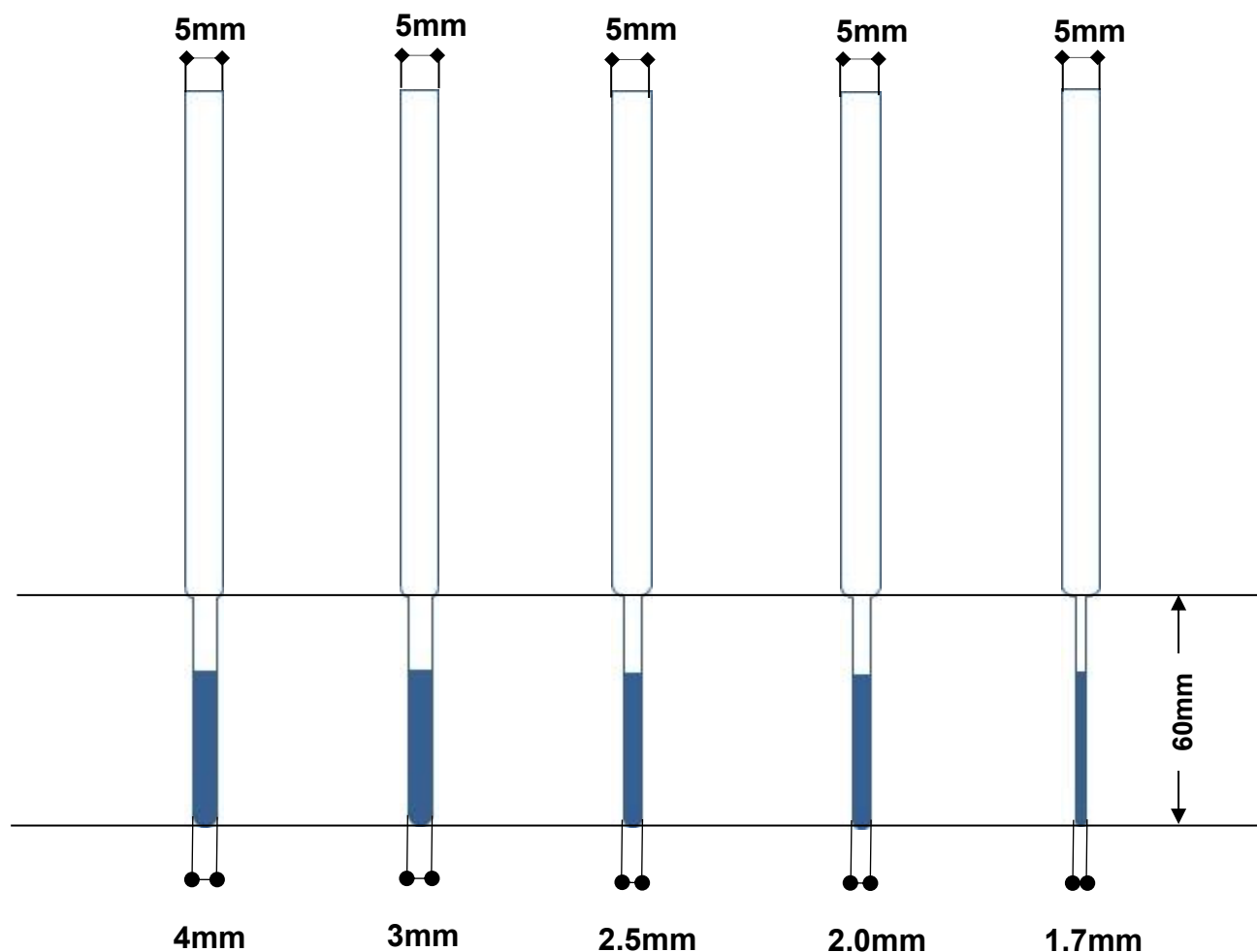

Volume comparison

Sample	4mm (SP-505)	3mm (SP-504)	2.5mm (SP-503)	2.0mm (SP-502)	1.7mm (SP-501)	5mm Thin wall
30mm	240uL	145 uL	84 uL	53 uL	34 uL	410uL
40mm	321 uL	194 uL	113 uL	70 uL	45 uL	549uL


Parts N.O.	Overall Length	Upper O.D.	Stem Length	Stem O.D.	Stem I.D.	List Price(JPY)
SP-501	180mm	5mm	60mm	1.7mm	1.2mm $\pm$ 0.01mm	4,100
SP-502	180mm	5mm	60mm	2mm	1.5mm $\pm$ 0.01mm	4,100
SP-503	180mm	5mm	60mm	2.5mm	1.9mm $\pm$ 0.01mm	2,900
SP-504	180mm	5mm	60mm	3mm	2.5mm $\pm$ 0.01mm	2,900
SP-505	180mm	5mm	60mm	4mm	3.2mm $\pm$ 0.01mm	2,900



5mm ϕ Coaxial NMR tube set

When you do not want to mix sample with solvent or to compare with the standard sample
We strong recommend the items.

We sale an outer tube(5mmO.D.), an inner tube with cap.

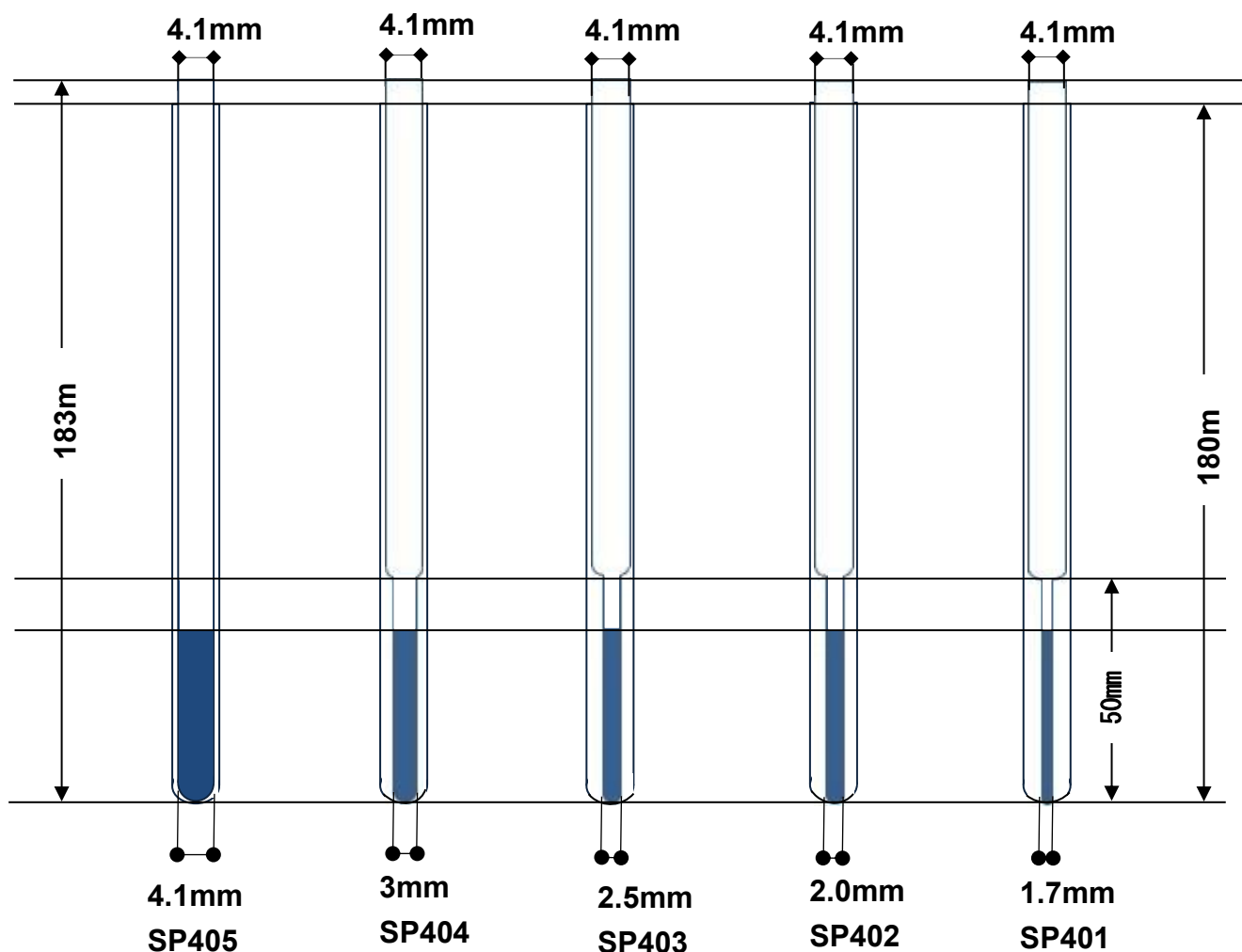

P.N.	Inner Tube Stem O.D.	Inner Tube Volume 40mm high	Outer Tube volume (When setting with an inner tube and an outer tube) 40mm
SP-401	1.7mm	45uL	459uL
SP-402	2mm	70uL	425uL
SP-403	2.5mm	113uL	355uL
SP-404	3mm	194uL	270uL
SP-405	4.1mm	317uL	55uL
SPT-401	1.7mm	45uL	547uL
SPT-402	2mm	70uL	512uL
SPT-403	2.5mm	113uL	442uL
SPT-404	3mm	194uL	257uL
SPT-405	4.1mm	317uL	142uL

5mmφ Coaxial NMR tube set

With PS-001-7(Thin wall Outer Tube O.D. 5mm)


Inner tube detail

P.N.	Length	O.D. Inner tube	Stem Length	Stem O.D	Stem I.D.	tolerance	1 Set(JPY)
SP-401	183mm	4.1mm	50mm	1.7mm	1.2mm	±0.01mm	6,800
SP-402	183mm	4.1mm	50mm	2mm	1.5mm	±0.01mm	6,800
SP-403	183mm	4.1mm	50mm	2.5mm	1.9mm	±0.01mm	5,600
SP-404	183mm	4.1mm	50mm	3mm	2.5mm	±0.01mm	5,600
SP-405	183mm	4.1mm	50mm	4.1mm	3.2mm	±0.01mm	5,600

With PST-001-7(Ultra thin wall Outer Tube O.D. 5mm)

Inner tube detail

P.N.	Length	O.D.	Stem Length	Stem O.D	Stem I.D.	tolerance	1 Set(JPY)
SPT-401	183mm	4.1mm	50mm	1.7mm	1.2mm	±0.01mm	7,900
SPT-402	183mm	4.1mm	50mm	2mm	1.5mm	±0.01mm	7,900
SPT-403	183mm	4.1mm	50mm	2.5mm	1.9mm	±0.01mm	6,800
SPT-404	183mm	4.1mm	50mm	3mm	2.5mm	±0.01mm	6,800
SPT-405	183mm	4.1mm	50mm	4.1mm	3.2mm	±0.01mm	6,800

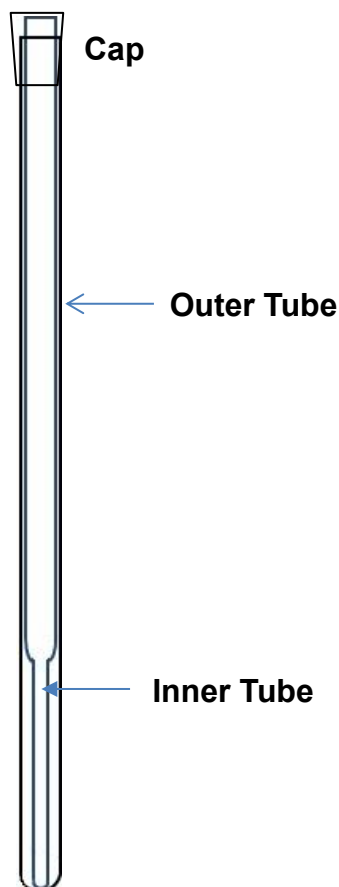

SPARE PARTS	JPY
SP-401INNER TUBE	4,100
SP-402INNER TUBE	4,100
SP-403INNER TUBE	2,900
SP-404INNER TUBE	2,900
SP-405INNER TUBE	2,900
PS-001 OUTER TUBE	2,200
PST-001 OUTER TUBE	3,500
CT-005(CAP)	700

We cannot guarantee for combinations with outer or inner tubes from other manufacturers.

5mm ϕ Coaxial INNER TUBE


for Heat sealing

The total length of the outer tube is 180mm, and the total length of this product is 200mm, so even if you fill the inner tube with sample or solvent and heat seal the top, it will still maintain a longer dimension than the outer tube.

We do not sell this as a set because it will not be possible to cap it. Please purchase the outer tube separately. PS-001-7, PST-001-7


☆We do not guarantee insertion into outer tubes from other manufacturers due to differences in inner diameter dimensions and tolerances.

P.N.	Length	Upper O.D.	Stem Length	Stem O.D.	Stem I.D.	tolerance	1 Peace (JPY)
SP-401L200	200mm	4.1mm	50mm	1.7mm	1.2mm	$\pm 0.01\text{mm}$	5,000
SP-402L200	200mm	4.1mm	50mm	2mm	1.5mm	$\pm 0.01\text{mm}$	5,000
SP-403L200	200mm	4.1mm	50mm	2.5mm	1.9mm	$\pm 0.01\text{mm}$	3,500
SP-404L200	200mm	4.1mm	50mm	3mm	2.5mm	$\pm 0.01\text{mm}$	3,500
SP-405L200	200mm	4.1mm	50mm	4.1mm	3.2mm	$\pm 0.01\text{mm}$	3,500

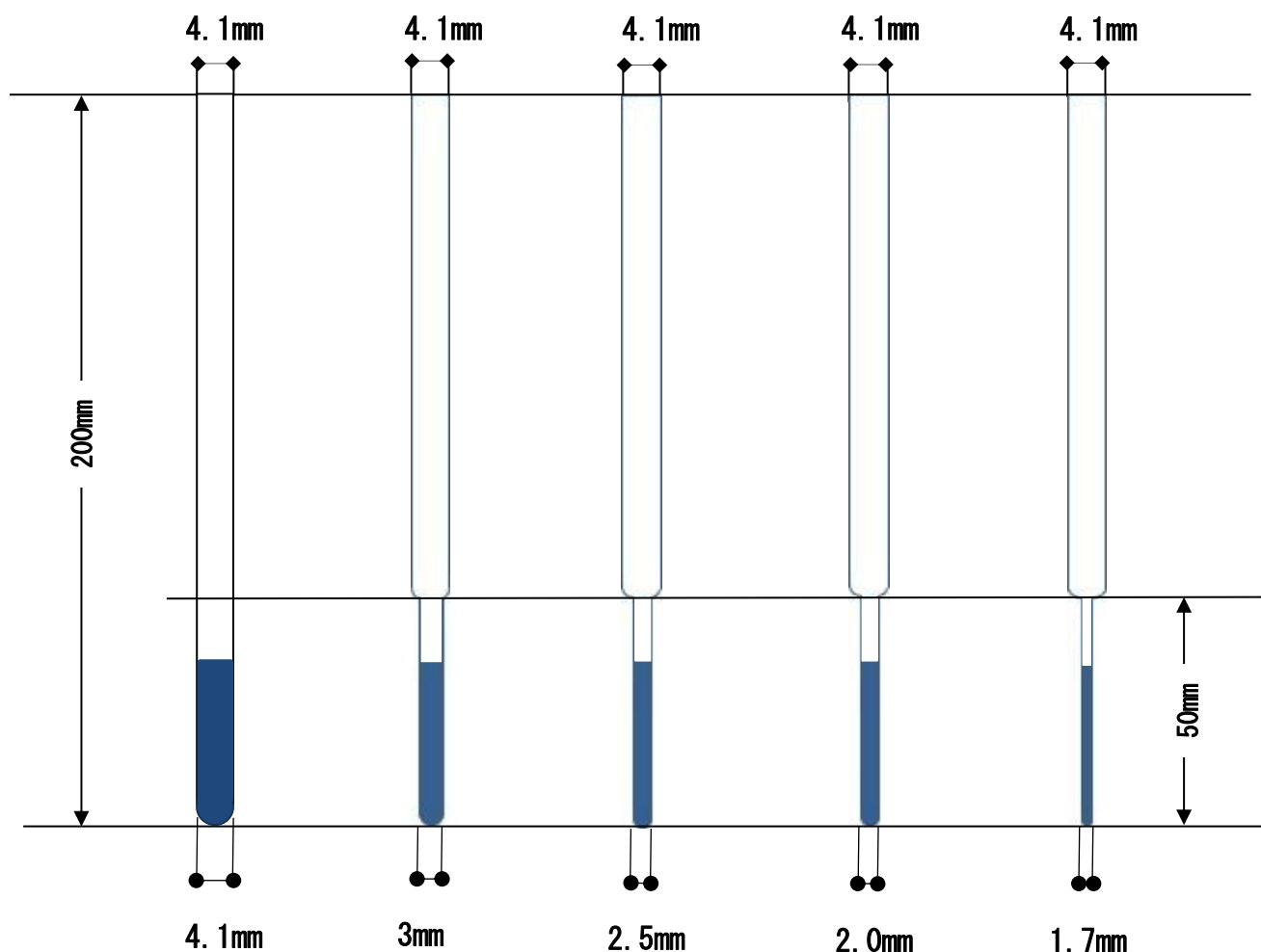

5mm ϕ Coaxial GAS tight NMR tube set

A coaxial tube that can block outside air. It is effective for samples that are sensitive to vacuum, nitrogen substitution, evaporation, water absorption, etc. Apply vacuum grease to the edge of the lid before use. Be careful not to let grease get into the sample tube. It can withstand long-term sealing.

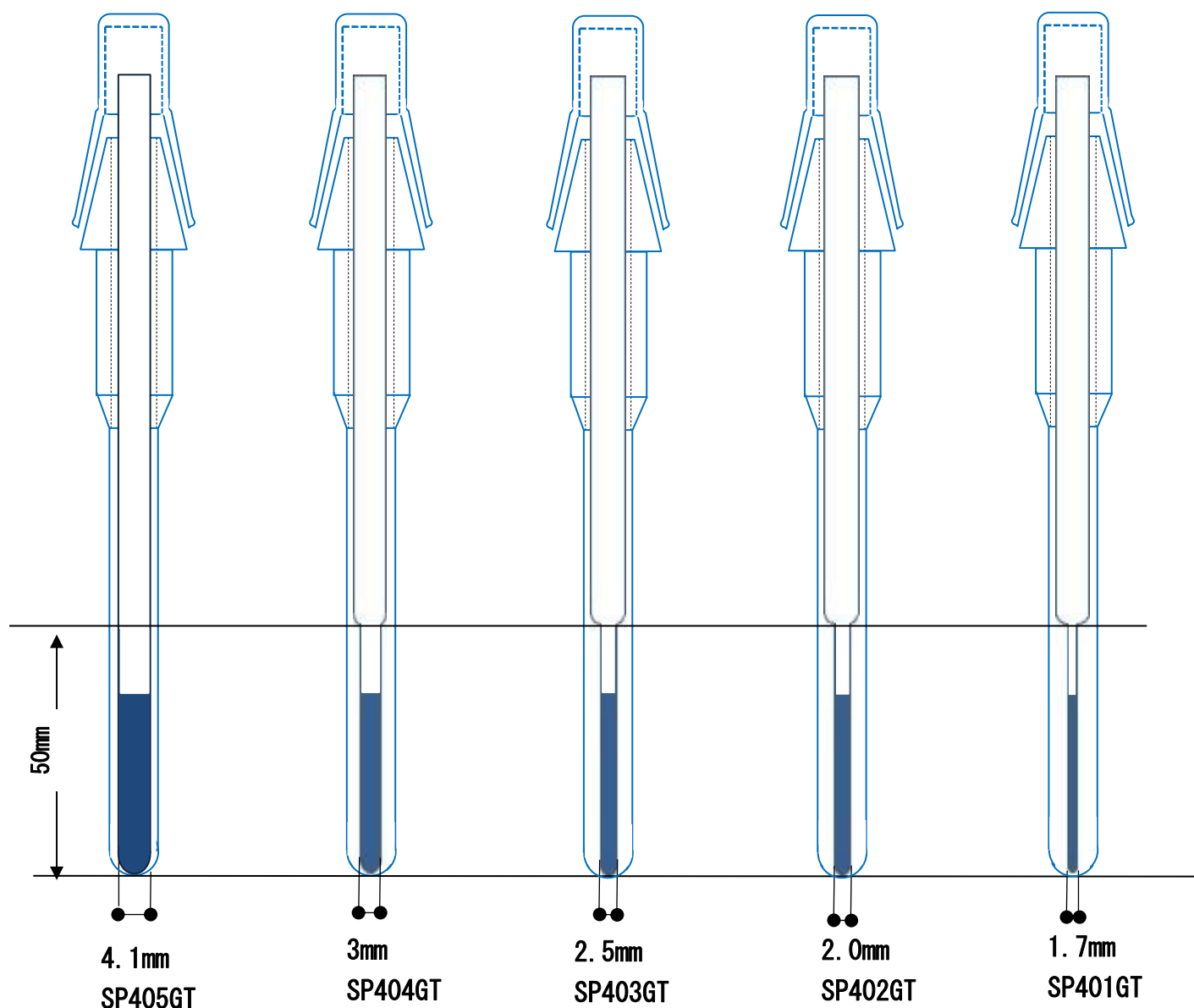

With PS-001-7GT(Thin wall GAS tight Outer Tube O.D. 5mm)

Inner tube detail

P.N.	O.D. Inner tube	Stem Length	Stem O.D	Stem I.D.	tolerance	1 Set(JPY)
SP-401GT	4.1mm	50mm	1.7mm	1.2mm	±0.01mm	15,000
SP-402GT	4.1mm	50mm	2mm	1.5mm	±0.01mm	15,000
SP-403GT	4.1mm	50mm	2.5mm	1.9mm	±0.01mm	13,800
SP-404GT	4.1mm	50mm	3mm	2.5mm	±0.01mm	13,800
SP-405GT	4.1mm	50mm	4.1mm	3.2mm	±0.01mm	13,800

With PST-001-7GT(Ultra Thin wall GAS tight Outer Tube O.D. 5mm)


Inner tube detail

P.N.	O.D. Inner tube	Stem Length	Stem O.D	Stem I.D.	tolerance	1 Set(JPY)
SPT-401GT	4.1mm	50mm	1.7mm	1.2mm	±0.01mm	16,100
SPT-402GT	4.1mm	50mm	2mm	1.5mm	±0.01mm	16,100
SPT-403GT	4.1mm	50mm	2.5mm	1.9mm	±0.01mm	15,000
SPT-404GT	4.1mm	50mm	3mm	2.5mm	±0.01mm	15,000
SPT-405GT	4.1mm	50mm	4.1mm	3.2mm	±0.01mm	15,000

SPARE PARTS		JPY
SP-401INNER TUBE		4,100
SP-402INNER TUBE		4,100
SP-403INNER TUBE		2,900
SP-404INNER TUBE		2,900
SP-405INNER TUBE		2,900
PS-001GT GAS Tight OUTER TUBE		11,000
PST-001GT GAS Tight OUTER TUBE		12,100



Quartz NMR Tubes

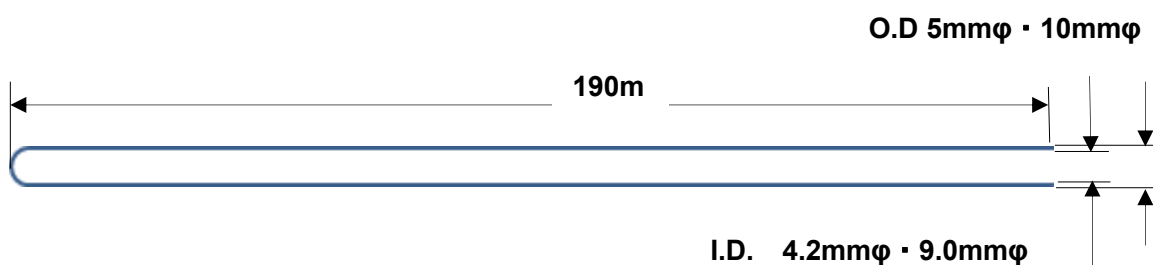

Dimension (mm)						
Parts N.O.	Length	O.D.	I.D.	Price/a peace	Material	For
SS-002	190	4.965 -0.005 -0.010	4.20 ±0.1	2,600	Fused Silica	NMR
SS-0010	190	9.98 ±0.01	9.0 ±0.1	4,600	Fused Silica	NMR


Fused Silica	Metallic impurity (ppm)										OH radical
	Na	K	Li	Mg	Mn	Fe	Ca	Al	Ti	Zr	
	1.3	<3.0	1.0	0.5	0.1	1.9	1.8	20.3	1.4	2.4	<5

Ultra thin wall NMR Tubes

Benefit


- ★ Tube thickness effect Signal to Noise
- ★ Volume gain
- ★ Sample approach RF coil
- ★ Better resolution of NMR spectra (especially Ultra High Field Magnet)
- ★ Easy to use

★ **25% gain Signal to Noise**

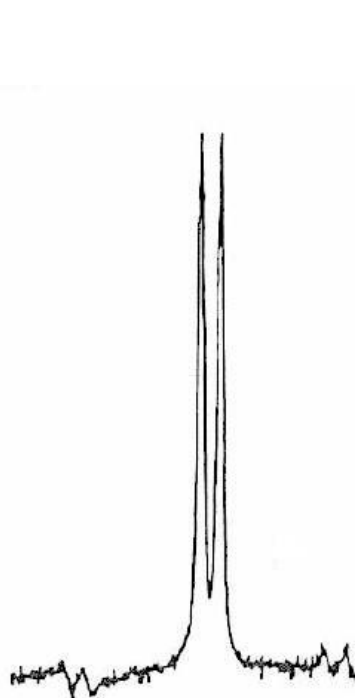

Ultra Precision NMR Sample Tubes

Figure 1

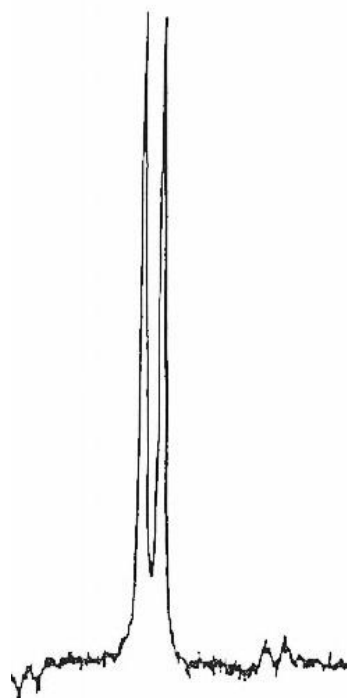

Thin wall NMR Tubes


Figure 2

¹H NMR Data 1 scan 20mM Sucrose in D₂O 298°K Bruker 600MHz

U.S. Patent No.5, 573, 567

Ultra Thin wall NMR Tubes



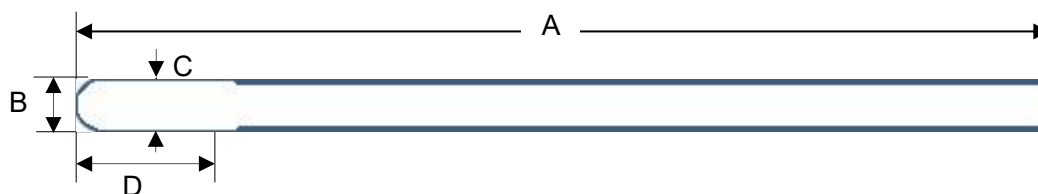


Parts N.O.	Dimension (mm)				Tolerance (μm)		Price/a peace (JPY)
	Length(A)	O.D.(B)	Wall Thickness(C)	Length of thin wall(D)	Wall thicknes s control	Straightnes s	
PST-001	180	4.965 +0.005 -0.010	0.21	50	10	10	3,500
PST-002	180	4.965 +0.005 -0.010	0.21	50	20	25	2,300
ST8-001	190	7.98 ±0.01	0.25	50	10	10	5,200
ST8-002	190	7.98 ±0.01	0.25	50	20	20	4,100
ST10-001	190	9.98 ±0.01	0.25	50	10	10	5,200
ST10-002	190	9.98 ±0.01	0.25	50	20	20	4,100

Volume comparison (40mm Height)

I.D(φ)	Thin wall	Ultra Thin Wall
5mm	550μL	637μL
8mm	1,507μL	1,703μL
10mm	2,465μL	1,712μL



Precision Large Volume NMR Tubes

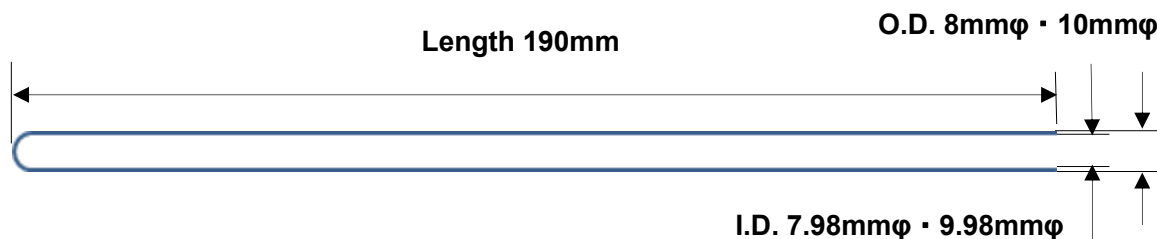

Parts N.O.	Dimension (mm)			Tolerance (μm)		Price/a peace (JPY)
	Length	O.D.	I.D.	Straightness	Wall thickness control	
S-8-001	190	7.98 ±0.01	7.03 +0.01 -0	10	15	3,500
S-8-002	190	7.98 ±0.01	7.03 +0.01 -0	10	20	2,900
S-10-001	190	9.98 ±0.01	9.03 +0.01 -0	10	15	3,500
S-10-002	190	9.98 ±0.01	9.03 +0.01 -0	10	20	2,900
S-10-003	190	9.98 ±0.01	9.03 +0.01 -0	20	30	2,300
S-10-004	190	9.98 ±0.01	9.03 +0.01 -0	20	40	2,100

If you need more volume, see “Ultra Thin wall (page 23)”


Volume comparison (40mm Height)

I.D(φ)	Thin wall	Ultra Thin Wall
5mm	550μL	637μL
8mm	1,507μL	1,703μL
10mm	2,465μL	1,712μL




ESR (EPR) Quartz Tubes

★Made of Fused Silica which is ultra-low background!!

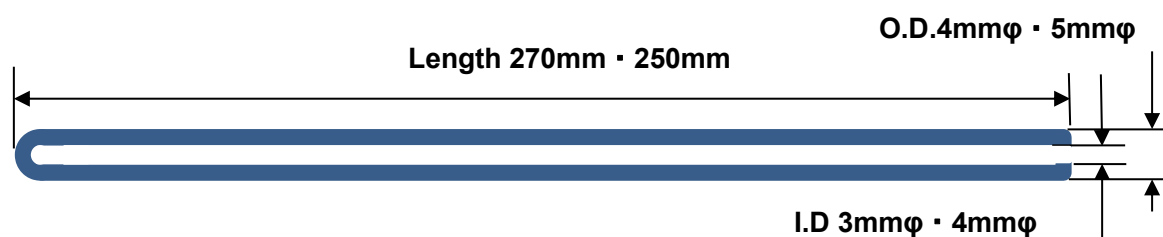

Part N.O.	O.D.	I.D.	Length	Material	Price/ a peace(JPY)
EPS-004J	4mm	3mm	270mm	Fused Silica	2,300
EPS-004B	4mm	3mm	250mm	Fused Silica	2,300
EPS-005J	5mm	4mm	270mm	Fused Silica	3,500
EPS-005B	5mm	4mm	250mm	Fused Silica	3,500

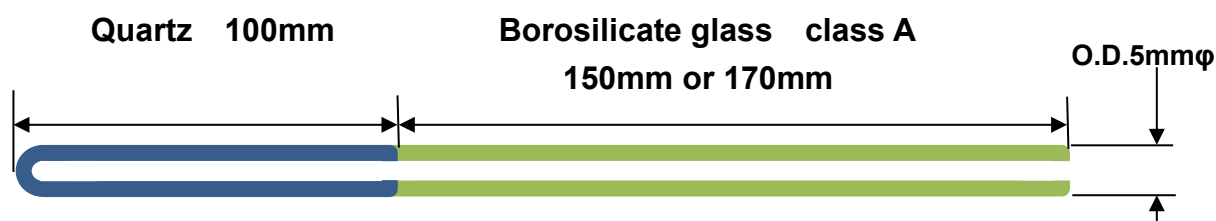

Part N.O.	O.D.	I.D(borosilicate glass)	I.D.(Quartz)	Length	Material	Price/ a peace(JPY)
EPW-005J	5mm	3mm	4mm	270mm	Quartz 100mm(Bottom) Borosilicate glass 170mm (Upper)	3,500
EPW-005B	5mm	3mm	4mm	250mm	Quartz 100mm(Bottom) Borosilicate glass 150mm (Upper)	3,500

The material produce by Shinetsu Quartz Co. LTD

	Metallic impurity (ppm)								OH radical
Fused Silica	Na	K	Li	Mg	Cu	Fe	Ca	Al	
HERALUX-E-LA	<0.05	<0.1	0.05	<0.05	<0.05	0.1	0.5	15	20


TRUE GAS TIGHT TUBE FOR NMR · ESR

It is effective for samples that are sensitive to vacuum, nitrogen substitution, evaporation, water absorption, etc. Apply vacuum grease to the edge of the lid before use. Be careful not to let grease get into the sample tube. It can withstand long-term sealing.

Benefit



1. To keep long term vacuum.
2. No need HEAT SEALING.
3. Standard Taper

§

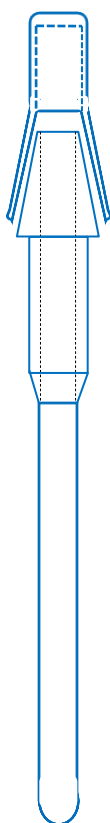

1. To keep long term vacuum.
2. No need HEAT SEALING.
3. Standard Taper

§

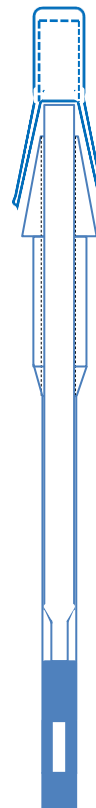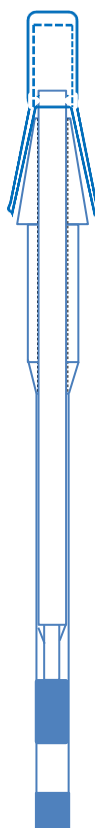

	P.N		JPY
5mm NMR Gas tight tube	PS-003GT	1Set	8,100
5mm symmetrical MICRO NMR Gas tight Tube for D ₂ O	BMS-005BGT	1Set	16,100
	BMS-005JGT	1Set	16,100
	BMS-005VGT	1Set	16,100
5mm symmetrical MICRO NMR Gas tight Tube for CDCl ₃	CMS-005BGT	1Set	16,100
	CMS-005JGT	1Set	16,100
	CMS-005VGT	1Set	16,100
5mm symmetrical MICRO NMR Gas tight Tube for DMSO((CD ₃) ₂ SO),Benzene (C ₆ D ₆), Pyridine (C ₅ D ₅ N)	DMS-005BGT	1Set	17,300
	DMS-005JGT	1Set	17,300
	DMS-005VGT	1Set	17,300
5mm symmetrical MICRO NMR Gas tight Tube for Methanol(CD ₃ OD),Acetonitrile (CD ₃ CN)用	MMS-005BGT	1Set	17,300
	MMS-005JGT	1Set	17,300
	MMS-005VGT	1Set	17,300
SLOT TUBE Gastight	SBMSGT	1Set	34,500
5mm NMR Gas tight tube(Quotz)	SS-002GT	1Set	10,400
5mm ESR Gastight tube (Quotz) length 250mm	EPS-005BGT	1Set	10,400
5mm ESR Gastight tube (Quotz) length 270mm	EPS-005JGT	1Set	9,200
4mm ESR Gastight tube (Quotz) length 250mm	EPS-004BGT	1Set	10,400
4mm ESR Gastight tube (Quotz) length 270mm	EPS-004JGT	1Set	9,200



Accessories

NMR Tube Caps



Parts N.O.	Description	Color	Price/100pieces
CP-005-Y	5mm	Yellow	1,800JPY
CP-005-G	5mm	Green	1,800JPY
CP-005-R	5mm	Red	1,800JPY
CP-005-P	5mm	Pink	1,800JPY

PTFE NMR Tube Caps

Parts N.O.	Description	Price/ a peace
CT-003	3mm	1,000JPY
CT-004	4mm	1,000JPY
CT-005	5mm	700JPY
CT-008	8mm	1,100JPY
CT-010	10mm	1,100JPY



FKM NMR Tube Caps

Parts N.O.	Description	Price/ a peace
CF-004	4.1mm	500JPY
CF-005	5mm	500JPY
CF-010	10mm	600JPY



FKM NMR Tube Caps

Parts N.O.	Description	Price/ a peace
CS-005	5mm	700JPY



5mm NMR SAMPLE TUBE for SampleJet

SJ-5

BENEFIT

- ☆High precision O.D. good for spinning, put on caps, insert spinner.
- ☆Material:DURAN ASTM ClassA good for variable temperature.
- ☆To melt end face for strong damage.



SPECIFICATION

PartNumber	(mm)				straightness(μm)
	Length	O.D.		I.D	
SJ-5	103.5(4inch)	4.965 +0.005 -0.010		4.220 ±0.10	30

Price

10packs(1,000PCS) (Minimum order unit)

10packsX1pack(100 pcs of tubes)(13,000JPY)=130,000JPY

CAPS for SJ-5

Just fit for SJ5 Price 1pack in 100 PCS of caps 5,800JPY



Part Number

SJC-5P(pink), SJC-5Y(yellow), SJC-5G(green), SJC-5W(white), SJC-5B(blue), SJC-5R(red)